

SpO2 Sensor Data Sheet

SPECIFICATIONS

> Infrared emitter*

- >Peak emission: 950nm
- >Centroid wave: 940nm
- >Spectral bandwidth: 42nm
- >Radiant intensity: 2mW/sr

> Red emitter*

- >Peak emission: 660nm
- >Centroid wave: 655nm
- >Spectral bandwidth: 17nm
- >Radiant intensity: 2.6mW/sr

> Detector*

- >Wavelength of max sensitivity: 920nm
- >Range of sensitivity: 400nm-1100nm
- >Radiant sensitive area: 1.3x1.3 (mm)
- >Spectral sensitivity(940nm): 0.77A/W

> Infrared/Red emitter

- >Duty cycle: 25%
- >Min current: 0.20mA
- >Max current: 50mA

> Resolution: 16bit

> Sampling frequency: 500Hz

FEATURES

- > Adjustable current for each LED using the API
- > Subtracts ambient light
- > Pre-conditioned digital output
- > High signal-to-noise ratio
- > Medical-grade raw data output
- > Ready-to-use form factor

APPLICATIONS

- > Oximetry
- > Heart rate & heart rate variability
- > Life sciences studies
- > Biomedical research
- > Human-Computer Interaction

GENERAL DESCRIPTION

The SpO2 (peripheral capillary oxygen saturation) sensor uses two emitting LED's one in the red region and the other in the infrared region of the spectrum. The reflected light of each one of these LED's is absorbed by a photodiode that converts this current into a digital value that is sent via SPI. This sensor can be used to estimate the oxygen saturation level



Fig. 1. Miniaturized form factor for minimally-intrusive application on the subjects

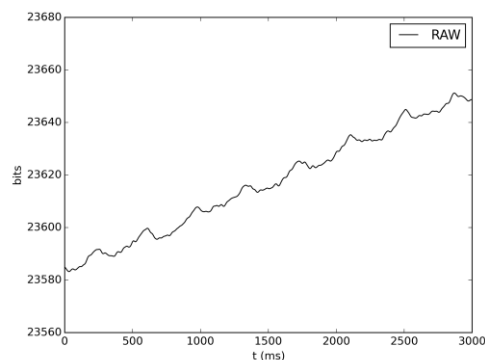


Fig. 2. Typical reflected red light raw data (acquired with biosignalsplux).

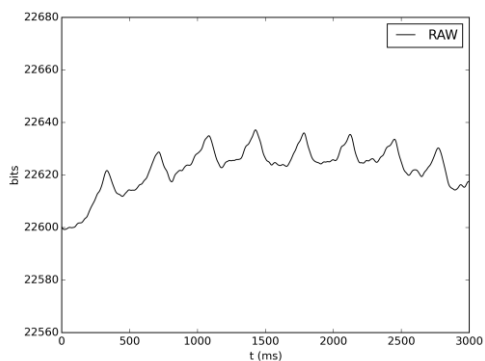


Fig. 3. Typical reflected infrared light raw data (acquired with biosignalsplux).

biosignalsplux
wearable body sensing platForm

PLUX – Wireless Biosignals, S.A.
Av. 5 de Outubro, n. 70 – 8.
1050-059 Lisbon, Portugal
plux@plux.info
<http://biosignalsplux.com/>

REV A

© 2015 PLUX

This information is provided "as is," and we make no express or implied warranties whatsoever with respect to functionality, operability, use, fitness for a particular purpose, or infringement of rights. We expressly disclaim any liability whatsoever for any direct, indirect, consequential, incidental or special damages, including, without limitation, lost revenues, lost profits, losses resulting from business interruption or loss of data, regardless of the form of action or legal theory under which the liability may be asserted, even if advised of the possibility of such damages.

* For additional parameters please contact plux@plux.info

SpO2 Sensor


Sensor Data Sheet

on the blood with +/- 2% accuracy compared to a medical sensor.

PHYSICAL CHARACTERISTICS


Infrared emitter relative spectral emission

$I_{rel}=f(\lambda)$, $T_A=25^\circ\text{C}$, $I_F=20\text{ mA}$

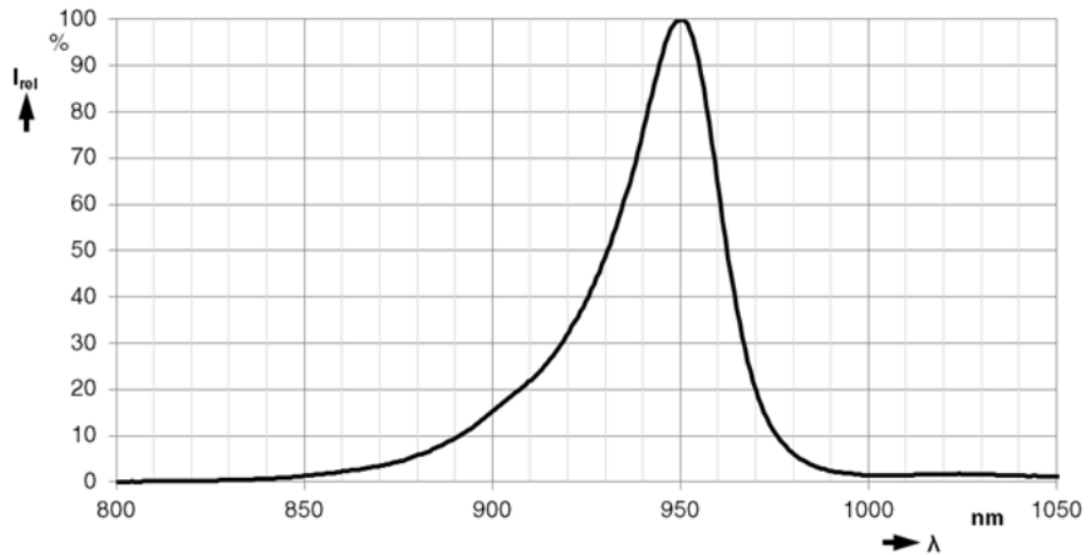

Red emitter relative spectral emission


$I_{rel}=f(\lambda)$, $T_A=25^\circ\text{C}$, $I_F=20\text{ mA}$

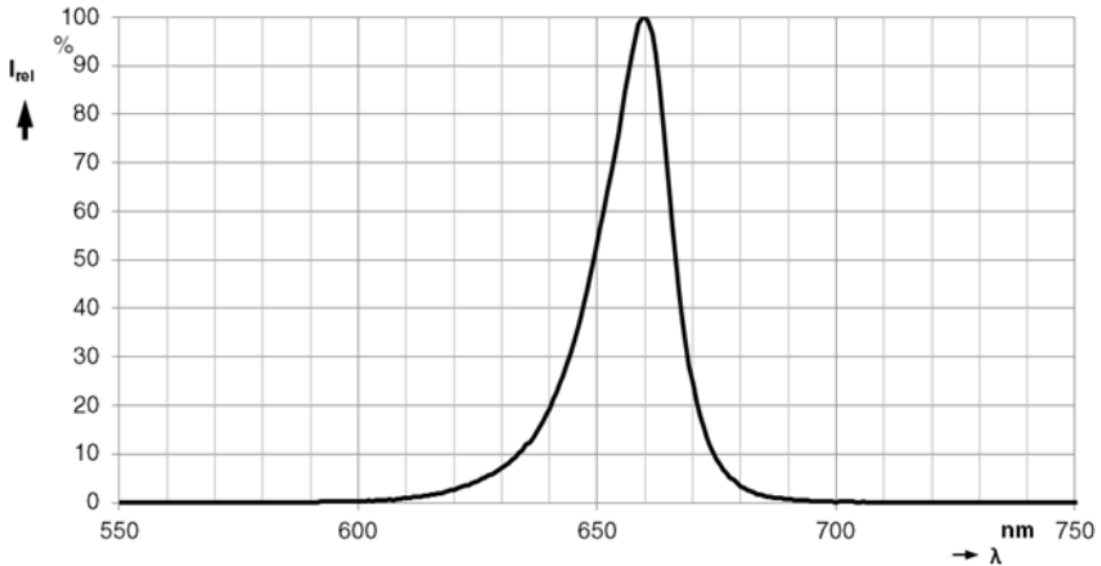

SpO2 Sensor Sensor Data Sheet

Detector relative spectral sensitivity

$S_{rel}=f(\lambda)$, $T_A=25^{\circ}\text{C}$

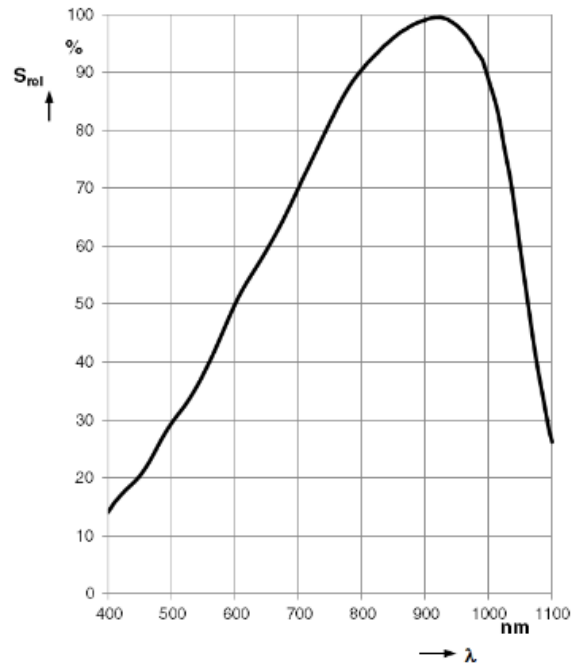

Function used to convert the photodiode current to a digital value:

$$V_{digital} = \frac{I_{pd} \times G}{1.2V \times 2^n}$$

I_{pd} : Photodiode current in Amperes
 G : Transimpedance Gain (1M Ω)
 n : Number of bits (8 or 16)

ORDERING GUIDE

Reference	Package Description